Cellpod MC:WELL

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Bert Sakmann Facts

Photo from the Nobel
Foundation archive.

Bert Sakmann
The Nobel Prize in Physiology or Medicine 1991

Born: 12 June 1942, Stuttgart, Germany

Affiliation at the time of the award: Max-Planck-Institut für
medizinische Forschung, Heidelberg, Federal Republic of
Germany

Prize motivation: "for their discoveries concerning the
function of single ion channels in cells."

Prize share: 1/2

Work

One of the fundamental processes of life is the transit of charged atoms - ions - through the surface layer of cells that make up organisms. Among other things, this transit is of crucial importance for the transfer of signals in nerves and muscles so that the body functions. Around 1980 Bert Sakmann and Erwin Neher developed a method for measuring the extremely weak currents involved in ion transits. These results confirmed that the transit occurs through ion channels - molecules on the surface of cells that under certain conditions allow ions to pass through.

PATCH CLAMP TECHNIQUES FOR STUDYING IONIC CHANNELS IN EXCITABLE MEMBRANES

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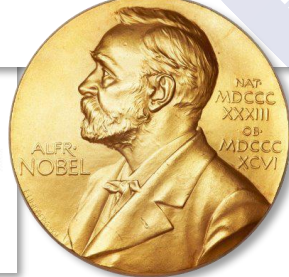
INTRODUCTION

Availability of techniques largely influences the choice of preparations in electrophysiological research. To study basic mechanisms of electrical excitability, usually the squid axon or giant snail neurons are used for technical reasons, or more simply, because they are large. For other reasons one would very often prefer to use mammalian cells, but they are much smaller and more difficult to handle. Recently, more freedom of choice was introduced by the so-called patch clamp techniques, which resulted from the technical challenge that the measurement of single-channel currents presented. It is evident that the smallness of these quantities required that the measurement be restricted to very small structures—small patches of membrane or very small cells. This led to the development of new techniques well suited for studying membrane currents from cells of any size and particularly suited to study small cells in culture. This review aims to describe the uses of the technique and to emphasize those results that use features specific to the various configurations of the patch clamp. We do not intend to present technical details, which can be found in the original literature (32, 74).

BASIC PRINCIPLE AND VARIANTS OF THE PATCH CLAMP METHOD

The original goal was to resolve the current contributions of individual ionic channels in a biological membrane. This was achieved by placing heat-polished

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